


Xmas Edition

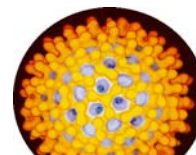

# APP-CHEM NEWSLETTER



**BEAT THE CREDIT CRUNCH!**

**EXTRA 10% DISCOUNT ON APP-CHEM REAGENTS UNTIL 22<sup>nd</sup> DECEMBER**

Simply quote "NL006" when ordering



## 2008 – A year of Investment and Re-growth

This Year App-Chem has seen continued growth in sales and this has enabled the management team to reinvest in the company; providing new laboratory equipment, fume cupboards, IT software and equipment and vocational staff training.

This investment has meant that our efficiency is even better than last year; with an average turnaround for reagents of just one day.

The AquaTech Exhibition in Amsterdam has already led to the appointment of three new distributors, with a further four lined up for the first quarter of 2009. Further announcements regarding these will be made in the New Year.

For 2009 we are re-launching the App-Chem brand with a new logo, new look catalogue and price lists and a **BRAND NEW** multifunctional website.

The website will facilitate the complete integration of the App-Chem system and will have:

- **Online Ordering**
- **Multi-lingual Safety Data Sheets**
- **Method Sheets and Instructions**
- **Technical Hints and Tips**
- **Regular Newsletters**

This will be launched early in the New Year.

## Rapid *Legionella* Field Test



The test kit contains all the items required to perform 5 tests – including filters, recovery buffer, collection vessel and sample point.

- These tests are optimised for use with **industrial** water systems, **cooling towers** and whirlpool **spas**.
- The filtration step needs **no pumping**, just connect the sample point to the water system and use the system pressure to drive the filtration.
- Single use **hollow fibre filtration modules**. Pre-filled syringe containing recovery buffer; no need to measure or dilute solutions.
- The test is **simple to use** and **fast**, with no interpretation required.
- Results are **easy to read** and obtained within **30 to 35 minutes**.
- The test allows **immediate understanding** and response to the water conditions.
- Use can **reduce risk** of outbreaks and litigation.
- *Legionella* control case



## Christmas Opening Hours

App-Chem will be closed from 1pm on 24<sup>th</sup> December 2008 and will not reopen until Monday 5<sup>th</sup> January 2009.

The staff at App-Chem would like to thank you for your continued support and look forward to dealing with you in 2009.



## Getting in touch

Whether you want to place an order, request a catalogue, have a query, or would like to arrange a visit please get in touch with us in one of the following ways.

### UK OFFICE

App-Chem House  
Hugmore Lane  
Llan y Pwll  
Wrexham  
LL13 9YE



### TELEPHONE/ FAX

TELEPHONE: +44 (0) 1978 661551

FAX: +44 (0) 1978 661611

graham.allen@app-chem.com  
catherine.bennion@app-chem.com  
aimee.williams@app-chem.com  
wendy.evans@app-chem.com  
susan.bentley@app-chem.com

Managing Director  
Technical Sales Manager  
Sales Administrator  
Sales Administrator  
Accounts Manager

[www.app-chem.com](http://www.app-chem.com)